

# GenePasQ Data Sheet / Reading guide

Title \_\_\_\_\_

DATE (                      )

Name (                      )

• Gene amplification by PCR


Multiplex PCR plus Kit

(Lot #) \_\_\_\_\_


①	95 °C	5 min	} 35 cycles
②	95 °C	15 sec	
③	60 °C	30 sec	
④	72 °C	15 sec	
⑤	4 °C		

• GenePasQ

(Lot #) \_\_\_\_\_

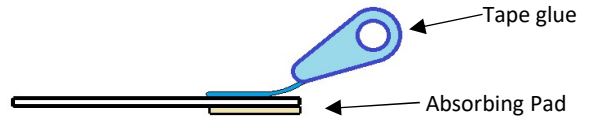

Paste the GenePasQ strips which is glued behind of absorbing pad with Tape glue.

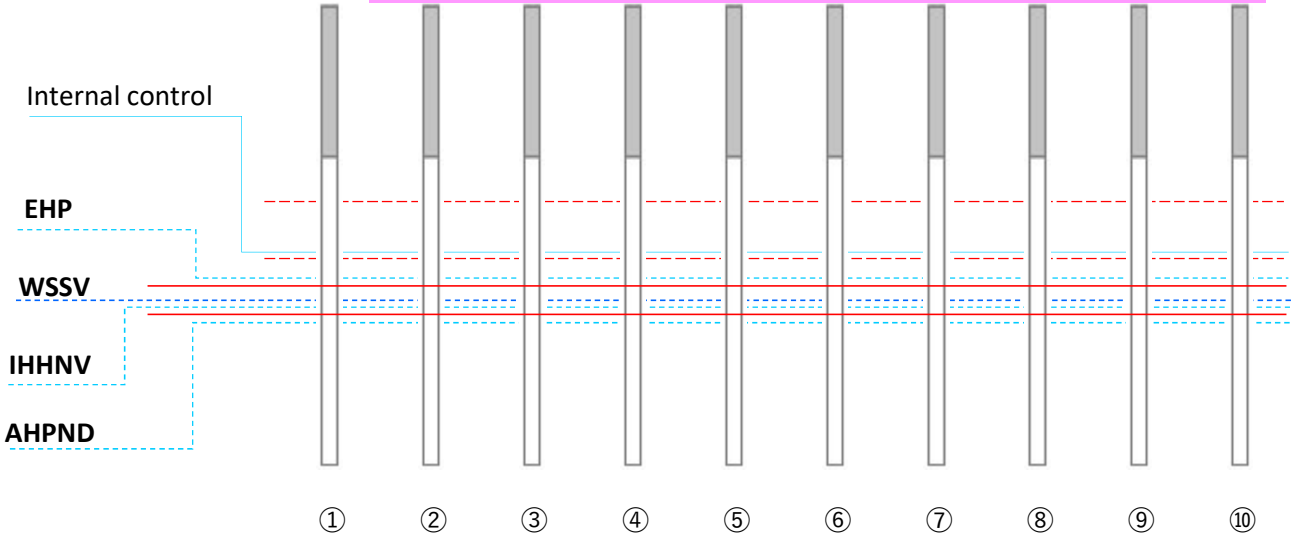

EHP	①	②	③	④	⑤	⑥	⑦	⑧	⑨	⑩
WSSV										
IHHNV										
AHPND										

\* Internal control (IC) line often becomes faint when a target gene is amplified.

test	comment
①	
②	
③	
④	
⑤	
⑥	
⑦	
⑧	
⑨	
⑩	